

Eichrom 10-50 mL Vial Shipping Systems Operation Manual

Part Numbers: SH-SYS-VIAL-30ML, SH-SYS-VIAL-30ML-LT, SH-SYS-VIAL-50ML-T

DIRECTIONS FOR USE

1. With the Vial Shield secure, twist the top counterclockwise to unscrew the top.
2. Place the vial with isotope inside the Vial Shield.
3. Ensure Vial Shield Rubber Seal is present and in good condition.
4. Carefully replace the Vial Shield top, twisting clockwise to tighten.
5. Place the Shield inside the cylindrical cavity of the Vial Shield Case.
6. Close the Vial Shield Case. To lock, place each of the latches onto its respective catch plate. Turn the latch handle to tighten. Flip latch handle down.
7. Seal Vial Shield Case with a zip tie or wire seal. Feed the wire or tie through spring and hole in the latch handle. Tighten wire or tie to prevent rotation of the latch handle.

NOTE: Seal may be optional when shipping in a sealed cargo compartment of a closed transport vehicle in exclusive use. Refer to applicable regulatory requirements.



Figure 1: Vial Shield Top, Vial Shield Bottom with Rubber Seal and Vial Inside, Vial Shield Case - Open



Figure 2: Vial Shield Case- Open with Vial Shield Inside



Figure 3: Vial Shield Case - Closed with Latches Locked. Zip Tie Seal Fed Through Spring and Hole in Latch Handle



Figure 4: Zip Tie Seal Tightened to Prevent Rotation of the Latch Handle